



CAR発現測定の ソリューションサービス

PE標識タンパク質 FITC標識タンパク質 Biotin標識タンパク質

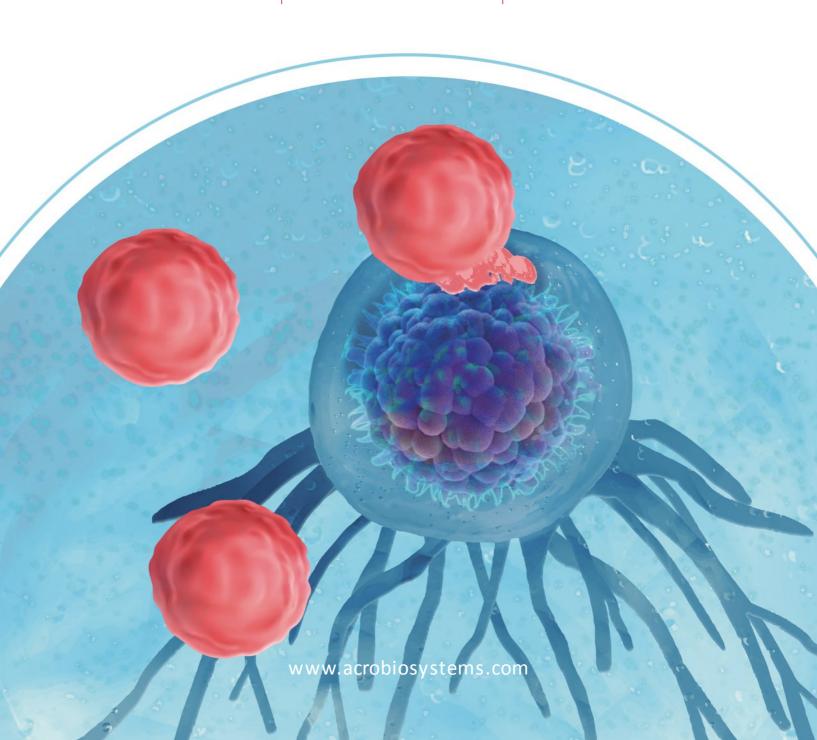

- フォーカスされている50近くのCAR-Tターゲット
- CAR発現の測定のために設計されている
- Protocolは無料で提供

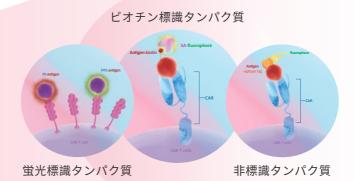





■ 確率したプラットホーム、 厳しい品質管理、品質管 理基準によって認定されている



Food and Drug Administration Silver Spring, MD 20993

DMF 034936

DMF ACKNOWLEDGEMENT

- ACROBIOSYSTEMS CO., LTD. FLOOR 4, BUILDING 4, No.8 HONGDA NORTH ROAD, BDA, BEIJING, CHINA
- 一部の製品は米国FDA DMFに登録済みです
- IND、NDAとBLAへの申請をサポートいたします
- DMF番号を直接使用できるので、製品の審査と 評価にかかる時間を節約できます


フォーカス製品



蛍光標識抗FMC63抗体によってAnti-CD19 CARの臨床サンプルを高感度で測定

製品コード	FM3-HPY53, FM3-FY45
製品タイプ	FITC、PEなどの様々な蛍光標識を提供しておる。
特長	FMC63 scFvの抗原認識部位を特異的に認識できる。特異性が強く、感度も高くて、臨床サンプルの測定に最適である。
お薦めのプラットホーム	ストリーミング分析プラットホーム
提供できるソリューションサービス	臨床サンプルには細胞の構成成分が複雑で、CAR-T細胞の含有量が低く、非特異的なバックグラウンドが強いといった問題があるため、これらの問題点を解決するサービスを提供しておる。

● PE標識抗FMC63抗体によってAnti-CD19 CARの臨床サンプルを高感度で測定

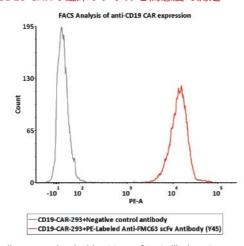


1e6 of the Anti-CD19 CAR-293 cells were stained with 100 µL of 1:50 dilution (2 µL stock solution in 100 µL FACS buffer) of PE-Labeled Monoclonal Anti-FMC63 scFv Antibody, Mouse IgG1 (Cat. No. FM3-HPY53) and negative control antibody respectively, PE signal was used to evaluate the binding activity.

蛍光標識CD19タンパク質によってAnti-CD19 CARの発現を特異的に測定

製品コード	CD9-HP2H3, CD9-HF2H2, CD9-HF251
製品タイプ	FITC、PEなどの蛍光標識を提供しておる。
特長	Anti-CD19 CAR-T細胞の品質確認に利用でき、Anti-CD19 CARの発現の測定も促進できる。
お薦めのプラットホーム	ストリーミング分析プラットホーム
提供できるソリューションサービス	非特異的なバックグラウンドが強い問題点を解決できるので、実験時間を半分に短縮できる。

➡ PE標識CD19によってAnti-CD19 CARの発現を特異的に測定

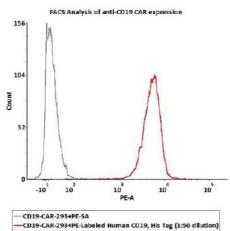


1e6 of the anti-CD19 CAR-293 cells were stained with 100 μ L of 1:50 dilution (2 μ L stock solution in 100 μ L FACS buffer) of PE-Labeled Human CD19 (20-291), His Tag (Cat. No. CD9-HP2H3). PE Streptavidin was used as negative control.




Star stainingは新世代の選択的蛍光標識タンパク質製品のシリーズである

製品コード	BCA-HF2H3、CD9-HF2H3
製品タイプ	FITC蛍光、BCMA、CD19などのターゲットも含まれている。
特長	新世代の選択的蛍光標識技術を用いることで、タンパク質の天然活性が保たれ、バッチ間の一貫性も高い。 CAR-T細胞の品質管理と臨床サンプル解析に設計されている。
お薦めのプラットホーム	ストリーミング分析プラットホーム
提供できるソリューションサービス	感度が低く、非特異的なバックグラウンドが強いなどの問題点を解決し、ワンステップで染色をし、実験時間を半分に短縮できる。

👈 Star Staining新世代の選択的蛍光標識BCMAタンパク質によってBCMA CARの発現を特異的に測定

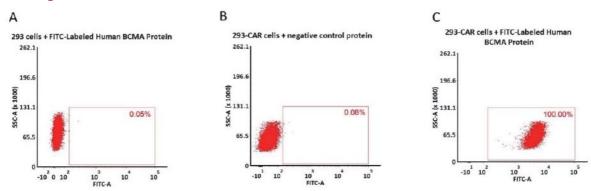

2e5 of 293 CAR cells transfected with anti-BCMA-scFv were stained with 100 μ L of 1 μ g/mL of FITC-Labeled Human BCMA, His Tag (Cat. No. BCA-HF2H3) and negative control protein respectively (Fig. C and B), and non-transfected 293 cells were used as a control (Fig. A), FITC signal was used to evaluate the binding activity.

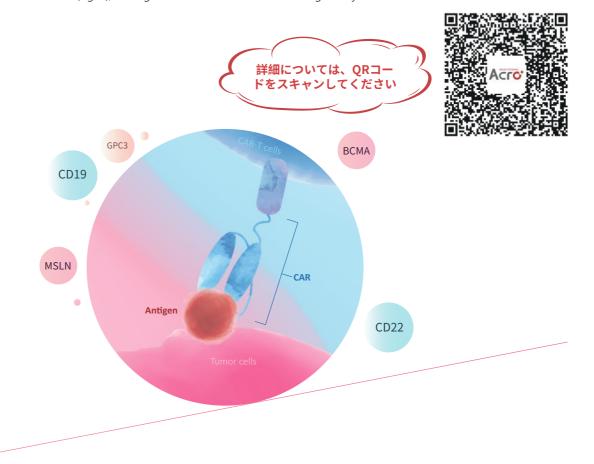

CAR-T標識タンパク質製品リスト(フォーカス製品)



♪ PE標識タンパク質製品リスト

分子名	製品コード	分子種	ラベル	発現システム
CD19	CD9-HP2H3	Human	His	HEK293
ВСМА	BCA-HP2H2 ^{NEW}	Human	His	HEK293
FMC63	FM3-HPY53 ^{NEW}	Human	Tag Free	HEK293
MSLN	MSN-HP2H5	Human	His	HEK293
SLAMF7	SL7-HP2H3	Human	His	HEK293
Siglec-3	CD3-HP2E3	Human	His	HEK293
CLEC12A	CLA-HP2Q3	Human	His	HEK293
Her2	HE2-HP2E3	Human	His	HEK293
CD30	CD0-HP2E3	Human	His	HEK293
CD4	CD4-HP2E3	Human	His	HEK293
CD7	CD7-HP2E3 ^{NEW}	Human	His	HEK293
EGF R	EGR-HP2E3	Human	His	HEK293
ROR1	RO1-HP2E3	Human	His	HEK293
Glypican 3	GP3-HP2E3 ^{NEW}	Human	His	HEK293
EGFRvIII	EGI-HP2E3	Human	His	HEK293
GUCY2C	GUC-HP2E3	Human	His	HEK293
CD5	CD5-HP2H6	Human	His	HEK293
PSMA	PSA-HP2Q3	Human	His	HEK293
FOLR1	FO1-HP2H9	Human	His	HEK293
B7-H3 (4Ig)	B7B-HP2H9	Human	His	HEK293

FITC標識タンパク質製品リスト

分子名	製品コード	分子種	ラベル	発現システム
BCMA	BCA-HF2H3 ^{Star staining}	Human	His	HEK293
	CD9-HF2H3 ^{Star staining}	Human	His	HEK293
CD19	CD9-HF2H2 ^{DMF}	Human	His	HEK293
	CD9-HF251 ^{DMF}	Human	Fc	HEK293
DCMA	BCA-HF254 ^{DMF}	Human	Fc	HEK293
BCMA	BCA-HF2H1 ^{DMF}	Human	His	HEK293
Her2	HE2-HF224	Human	His	HEK293
FMC63	FM3-FY45 ^{NEW}	Human	Tag Free	HEK293
CD22	SI2-HF2H6	Human	His	HEK293
GPC3	GP3-HF2H1	Human	His	HEK293
MSLN	MSN-HF223	Human	His	HEK293
EGF R	EGR-HF2H5	Human	His	HEK293
PSMA	PSA-HF244	Human	His	HEK293
GUCY2C	GUC-HF2H8	Human	His	HEK293
CD70	CDD-HF2D4	Human	His	HEK293

分子名	製品コード	分子種	ラベル	発現システム
CLEC12A	CLA-HF247	Human	His	HEK293
FAP	FAP-HF243 ^{NEW}	Human	His	HEK293
CD138	SD1-HF228	Human	His	HEK293
CD20	RIB-FY35c	Mouse	Tag Free	HEK293
B7-H3	B7B-HF2E7	Human	His	HEK293

• ほかのFITC標識タンパク質

CD38	Her3	CD4	ROR1	CEACAM5
CD33	CD30	CD123	IL13RA2	CD147
Nectin-4	VEGF R2	EpCAM	CD8	CD7
FOLR1	EGFRvIII	SLAMF7	CD5	NKG2D
CD99	CD98			

Biotin標識タンパク質製品リスト

分子名	製品コード	分子種	ラベル	発現システム
DCMA	BC7-H82F0 ^{DMF}	Human	Fc & Avi	HEK293
BCMA	BCA-H82E4 ^{DMF}	Human	His & Avi	HEK293
CD10	CD9-H82F7	Human	Fc & Avi	HEK293
CD19	CD9-H82E9	Human	His & Avi	HEK293
CD22	SI2-H82F8	Human	Fc & Avi	HEK293
FMC63	FM3-BY45	Mouse	Tag Free	Hybridoma
GPC3	GP3-H82E5	Human	His & Avi	HEK293
Her2	HE2-H82E2	Human	His & Avi	HEK293
	MSN-H82E9	Human	His & Avi	HEK293
MSLN	MSN-H82F6	Human	Fc & Avi	HEK293
CEACAM5	CE5-H82E0	Human	His & Avi	HEK293
EGF R	EGR-H82E3	Human	His & Avi	HEK293
	GUC-H82E9	Human	His & Avi	HEK293
GUCY2C	GUC-H82F8	Human	Fc & Avi	HEK293
HGF R	MET-H82E1	Human	His & Avi	HEK293
60100	SY1-H82E2	Human	His & Avi	HEK293
CD138	SY1-H82F3	Human	Fc & Avi	HEK293
	FAP-H82Q6 ^{NEW}	Human	His & Avi	HEK293
FAP	FAP-C82Q5 ^{NEW}	Cynomolgus	His & Avi	HEK293
	FAP-M82Q5 ^{NEW}	Mouse	His & Avi	HEK293
CA125	CA5-H82F4	Human	Fc & Avi	HEK293
CD70	TN7-H82F4	Human	Fc & Avi	HEK293
CLEC12A	CLA-H82E6	Human	His & Avi	HEK293
CLECIZA	CLA-H82F7	Human	Fc & Avi	HEK293

• ほかのBiotin標識タンパク質

CD38	Her3	Nectin-4	CD133	EGFRvIII
CD4	ROR1	CD147	EGFRvIII	EpCAM
FOLR1	SLAMF7	CD8	CD30	VEGF R2
CD33				

4

非標識タンパク質製品リスト

分子名	製品コード	分子種	ラベル	発現システム
CD10	CD9-H52H2 ^{DMF}	Human	His	HEK293
CD19	CD9-H5251 ^{DMF}	Human	Fc	HEK293
BCMA	BC7-H5254 ^{DMF}	Human	Fc	HEK293
BCMA	BCA-H522y ^{DMF}	Human	His	HEK293
FMC63	FM3-Y45	Mouse	Tag Free	Hybridoma
FIVICOS	FM3-Y45P1 ^{NEW}	Mouse	Tag Free	HEK293
GPC3	GP3-H5258	Human	Fc	HEK293
GPC3	GP3-H52H4	Human	His	HEK293
MSLN	MSN-H5253	Human	Fc	HEK293
MSLN	MSN-H522a	Human	His	HEK293
Her2	HE2-H5225	Human	His	HEK293
CD22	CD2-H52H8	Human	His	HEK293
EGF R	EGR-H5252	Human	Fc	HEK293
CEACAM-5	CE5-H5226	Human	His	HEK293
PSMA	PSA-H52H3	Human	His	HEK293
HGF R	MET-H5227	Human	His	HEK293
GUCY2C	GUC-H52H5	Human	His	HEK293
CLEC12A	CLA-H5266	Human	Fc	HEK293
FAP	FAP-H5244 ^{NEW}	Human	His	HEK293
CD138	SY1-H5225	Human	His	HEK293

• ほかの非標識タンパク質

CD38	CD33	CD70	IL-3 R alpha	CD8 alpha
Her3	Nectin-4	CD30	EpCAM	IL-13 R alpha 2
ROR1	PSCA	EGFRvIII	SLAMF7	NKG2D
FOLR1				

BIOSYSTEMS

Her2 BAFFR LAG-3 Fc Receptor Siglec-10 **Biotinylated Protein** PD-L1 VEGF165 CD3 epsilon PD-1BCMA CD27PVRIG CD47 PSMA **OFGL1TFPI**Siglec-15 Integrin CD24 CD3E & CD3D CD20 CD19 FcRn PCSK9 IL-2 R alpha **CAR-T Target Protein** Glypican 3Integrin 5 TOTAL Service TOTAL SER **Biotinylated Protein** 20CD200GITR & Nectin-4 VEGF165 CD69 FGLI VEGF165 PCSK9 IgG1 Fc CD40 ^{ਡੂ}PD-L1 SIRP alpha ADA Service SPR /BLI analytical service



+86 400-682-2521 (Asia & Pacific) +1 800-810-0816 (US/Canada) order@acrobiosystems.com www.acrobiosystems.com 1 Innovation Way, Newark, DE 19711



